

ARCS PROCEDURE: Author: A. Koontz	DEACTIVATING THE WSI FROM ADAM	PRO(WSI)-011.000 May 4, 1999 Page 1 of 1
--	-----------------------------------	--

Deactivating the WSI from ADaM

I. Purpose:

Describes the procedure for deactivating the WSI from ADaM.

II. Cautions and Hazards:

None.

III. Requirements:

None.

IV. Procedure:

A. Steps:

1. On the X-terminal press the Instruments icon.
2. Select the WSI instrument icon.
3. Select "Deactivate" from the command list.
4. Select "Execute command."
5. Click the mouse in the black area of the password block. The cursor should be blinking.
6. Enter the password **ADaM**.
7. Press OK.
 - If the password block times out before you can type the password, then you can restart at step 3.

V. References:

None.

VI. Attachments:

None.