

ARCS PROCEDURE:  Author: A. Koontz	DEACTIVATING AN INSTRUMENT FROM ADAM	PRO(ADAM)-018.000  May 6, 1999 Page 1 of 1
--	---	---

## Deactivating an Instrument from ADaM

### I. Purpose:

Describes the process for deactivating an instrument from ADaM.

### II. Cautions and Hazards:

None.

### III. Requirements:

None.

### IV. Procedure:

#### A. Steps:

1. On the X-terminal press the Instruments icon.
2. Select the desired instrument icon.
3. Select 'Deactivate' from the command list.
4. Select 'Execute command.'
5. Click the mouse in the black area of the password block. The cursor should be blinking.
6. Enter the password **ADaM**.
7. Press OK.
  - If the password block times out before you can type the password, then you can restart at step 3.

### V. References:

None.

### VI. Attachments:

None.