

ARCS PROCEDURE: Author: A. Koontz	ACTIVATING AN INSTRUMENT FROM ADAM	PRO(ADAM)-019.000 May 6, 1999 Page 1 of 1
--	---------------------------------------	---

Activating an Instrument from ADaM

I. Purpose:

To describe the process for activating an instrument from ADaM.

II. Cautions and Hazards:

None.

III. Requirements:

None.

IV. Procedure:

A. Steps:

1. On the X-terminal press the Instruments icon.
2. Select the desire instrument icon.
3. Select 'Activate' from the command list.
4. Select 'Execute command.'
5. Click the mouse in the black area of the password block. The cursor should be blinking.
6. Enter the password **ADaM**.
7. Press OK.
 - If the password block times out before you can type the password, then you can restart at step 3.

V. References:

None.

VI. Attachments:

None.